Contents:

- 1 Aluminium container with 100 test sticks
- 2 Bottles Arsenic-1
- 2 Bottles Arsenic-2
- 1 Reaction vessel
- 1 syringe 10 ml
- 1 needle

Hazards informations:

Each test field on the test stick contains mercury (II) bromide. Toxic by contact with skin.

Method of application:

1. Use syringe to add 2 x 10 ml sample solution into the reaction vessel .
2. Add 1 measuring spoon reagent Arsenic-1 and swirl reaction vessel gently.
3. Add 1 measuring spoon reagent Arsenic-2.
4. Remove only as many test sticks as are required, and reseal the container immediately after use. **Do not touch the test paper zone.** Insert test strip with test field 2 cm into reaction vessel and clamp it with lid (see sketch). For pressure balance pricking the needle through the lid and leave it there.
5. During 30 minutes reaction time swirl 3 x gently the reaction vessel. The test field should not get in contact with the sample.
6. After 30 minutes remove test strip from reaction vessel and dip it for 2 sec. with the test field into dist. water.
7. Shake off excess liquid and compare test field to colour scale.

Interferences:

The following ions only interfere in larger concentrations:
 1000 mg/l: Ca²⁺, Fe²⁺, K⁺, Mg²⁺, Na⁺, Zn²⁺, Cl⁻, CN⁻, CO₃²⁻, SO₄²⁻, EDTA.
 500 mg/l: Fe³⁺, F⁻; 100 mg/l: Al³⁺, Sn²⁺, CrO₄²⁻, NO₂⁻, NO₃⁻, PO₄³⁻; 25 mg/l: ClO₃⁻;
 10 mg/l: Ni²⁺; 5 mg/l: Co²⁺; 1 mg/l: Ag⁺, Sb³⁺, SeO₃²⁻, SO₃²⁻, S₂O₃²⁻; 0.5 mg/l: Cu²⁺, S²⁻.

Storage:

Avoid exposing the sticks to sunlight and moisture. store the kit below +30° C in a dry place.

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